

# Customer Support

# Software Release Bulletin

**Date of Issue: 03/10/2024**

**Product: AARI**

**Version: 1.7.4.0**

**Minor Release**

# AARI

## Version 1.7.4.0

### Release Type: **Minor**

This is a Minor software release to address some user requests alongside a number of bug fixes.

### Key Features

- Arabic Reporting
- Preset management tweaks
- Licence corrections (DCS6 compatibility)

### **IMPORTANT**

This version of software will require a new license generated by f+f to enable a successful install. Please email [CustomersupportTeam@fosterfreeman.com](mailto:CustomersupportTeam@fosterfreeman.com) requesting a new license. Please ensure you include the serial number of the device within the request and would recommend requesting the new license prior to arranging the upgrade, so you have it available. The software will

## Installation Instructions

1. Log into the Distributor Website or Customer Portal as appropriate, navigate to the relevant product area, select the Software Updates drop down and Click the 'Download' button. Screenshot for illustration purposes only.

# AARI®

Streamlined visualisation and mark-up of fingerprints: using AARI, tasks that previously took many hours of manual processing can now be achieved in minutes. [View product at fosterfreeman.com](http://fosterfreeman.com)

User Manuals	+
Software Updates	-


Please download and read the Software Release Bulletin before installing or updating the software on your device.

AARI Software Release Bulletin Version 1.5.3.13 (21-12-23) [DOWNLOAD](#)  
1 file(s) 326.12 KB

AARI SW Release v1.5.3.13 20-12-23 [DOWNLOAD](#)  
1 file(s) 451.07 MB

2. Right click the file and select "Properties"
3. At the bottom right of the property's windows tick the box that says "Unblock"
4. Click Apply, then Click OK
5. Right click the file and select "Extract All.."
6. Ensure "Show Extracted Files" is checked and click "Extract"
7. Where products are not on a desktop computer, copy the file onto the USB drive and transfer the file to a folder onto the target hardware
8. On the target hardware, open the containing folder and double-click the executable file
9. Follow the on-screen instructions
10. Once completed, restart the device

As part of the installation, you will be asked for a new license with a similar screen to the below.

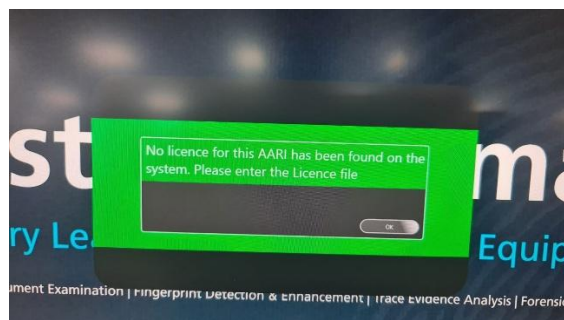

## Bug Fix / Feature Breakdown

ID	Title	Work Item Type
18331	Bug in new White Balance options on AARI	Change Request
20033	AARI Report not generated in Arabic	Change Request
20822	Ability to edit default presets	Change Request
20823	Ability to re-order image gallery by recent or oldest images	Change Request
28655	AARI Export image with all marked ROIs in independent file	Change Request
32121	AARI - Remember settings after re-opening software	Change Request
32122	AARI - progress indicator during export	Change Request
34337	AARI - License file incompatible with DCS6 license	Change Request
20538	Serial Number Improvements	Software Task
23146	ROI Region Original Image	Software Task
34657	'Delete preset ' button is available for default preset	Bug
34658	Overwriting presets is setting incorrect ISO values to the presets	Bug
34659	'Reset Preset' option is not available for some of the default presets	Bug