

Customer Support

Software Release Bulletin

Date of Issue: 21/08/2025

Product: Crimelite-AUTO / DISCOVER
Version no: 2.1.0.0

Minor RELEASE

CL-AUTO / DISCOVER

Version 2.1.0.0

Release Type: **Minor**

This is a Minor software release addressing numerous identified bugs and improvements as detailed in the table below.

Additional Feature Highlights

- Traditional Chinese Improvements
- Feature Improvement for DISCOVER
 - Exposure Control
 - Text Annotation

Software Download Instructions

Log into the Distributor Website or Customer Portal as appropriate, navigate to the relevant product area, select the Software Updates drop down and Click the 'Download' button next to the 'SW Release' to download the zip file. Below image for illustration purposes only.

Crime-lite[®] AUTO

Forensic imaging technology meets powerful multi-spectral illumination in this complete solution to the search, detection, and capture of evidence.

[View product at fosterfreeman.com](#)

User Manuals	+
Software Updates	-

Please download and read the Software Release Bulletin before installing or updating the software on your device.

CL AUTO & DISCOVER Software Release Bulletin v1.3.0 06-08-24 1 file(s) 340.25 KB	DOWNLOAD
CL-AUTO/DISCOVER SW Release v1.3.0 06-08-24 1 file(s) 533.71 MB	DOWNLOAD

Video Guides	+
--------------	---


Installation Instructions

This version can be installed directly over a previous version.


Please ensure you use a good quality NTFS formatted USB 3.0 stick on which to load the software update.

All CL-AUTO & DISCOVER software is installed via updates to the CL-AUTO unit itself.

1. Download the latest software via download link
2. Open zip file and extract AFU file. IE AUTO_2_0_0_0.AFU
3. Copy that AFU file onto a formatted (**NTFS**) USB drive. Make sure it is not inside any directory and also make sure it's the only update file.
4. Power up CL-Auto unit
5. Insert the USB 3.0 stick into the camera
6. Wait until you have access to power menu

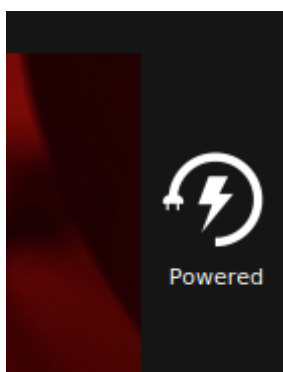

7. Open Settings menu

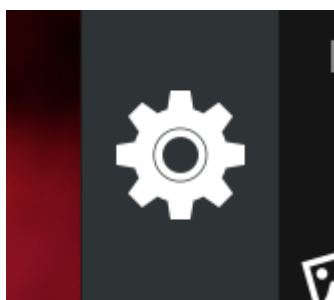

8. Click the "System and Update" option
9. Click "Install Update"
10. You should now see an install progress bar
11. Upon successful installation, the close button should turn into "Restart"
12. Select Restart to complete the installation

Some additional User Guide information related to a number of the new features has been provided at the end of this release note and is also available as a separate download from the portals.

This version contains the following enhancements and fixes.

ID	Change	Work Item Type
22362	DISCOVER CSU issues with zoom	Change Request
24428	Double click preview on DISCOVER to take a preview	Change Request
25834	When coax is inserted, OLM is activated automatically instead of coax on CL AUTO	Change Request
25911	Add Text annotation feature to DISCOVER	Change Request
30050	DISCOVER - Make 'Preview' button easier to select	Change Request
30051	DISCOVER - Add exposure compensation	Change Request
30833	Auto - not saving RAW after deleting images off SD card	Change Request
31428	Mode in Auto to export image only, without Hash/Metadata file	Change Request
28071	Exposure compensation slider replicated on Discover monitor?	Minor Project Task
32201	Traditional Chinese Language Review	Minor Project Task
32262	Exposure compensation value displayed on the Discover monitor	Minor Project Task
32266	Confirmation that SD card is safe to remove	Minor Project Task
32053	Add a TFT Pixel Test to the engineering screen	Software Task
14299	Discover - No difference between High and Low contrast presets if manual camera settings not selected	Bug
16962	Discover - Lifts preset - lighting options buttons text cut off	Bug
17534	Some light source data cut off on left hand menu using PVL presets	Bug
17545	Images captured with the App have blank thumbnails when viewed in App	Bug
17567	Discard button text in Notes is clipping in French, Russian, Spanish, German, Dutch, Italian	Bug
17673	Discover - Lifts preset text clipping in Polish, French and Spanish	Bug
22590	Coax and oblique light source data not shown in exported image PDF	Bug
22665	Discover images lost if monitor switched off/disconnected	Bug
23781	Unable to find RAW photos in SD Card	Bug
24075	Discover - Frequent USB issues	Bug
25071	Gallery thumbnails not visible in CL mobile app	Bug
31469	Discover - Mouse Disappears on re-insertion of USB dongle	Bug
32112	Measurement Tool - Zero allow as a measurement	Bug
32261	Exposure compensation not replicated across screens	Bug
32265	Option for exporting meta data not working	Bug
32323	Discover - Images in viewer names and positions change after switching monitor off then on	Bug
32325	Discover - Cursor jumps to bottom right corner of screen when any Preset option clicked	Bug
32326	Discover - language selection list items not being selected	Bug
32332	Traditional Chinese language cut off at bottom of screen	Bug
32338	Some images on SD cards are not showing metadata	Bug
32424	Image order changed in gallery after re-inserting SD card	Bug
32431	First time opening gallery after capturing an image not working	Bug

ID	Change	Work Item Type
32432	Secondary menu's overlapping primary menu	Bug
32578	Crash when coax USB connected	Bug
32579	Crash when oblique USB connected	Bug
32601	Discover - software sometimes logs out while adjusting intensity of peripheral light sources	Bug
32666	Custom Presets crashing/missing add button	Bug